

for
10mL
5mL
pipette

**No electric power
required at all!**

Repetitive Pipette

Peggy




Something Different.


WATSON BIO LAB
MADE IN JAPAN SINCE 1980



Aspirate the sample all at once, dispensing its necessary volume repetitively.

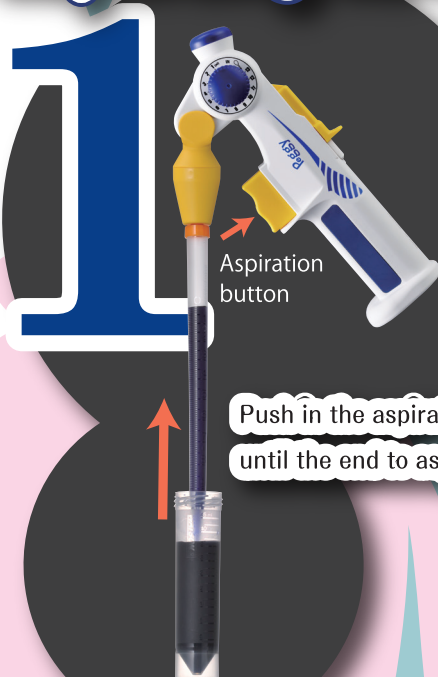

Push in the aspiration button once slowly
until the end to aspirate the set volume.

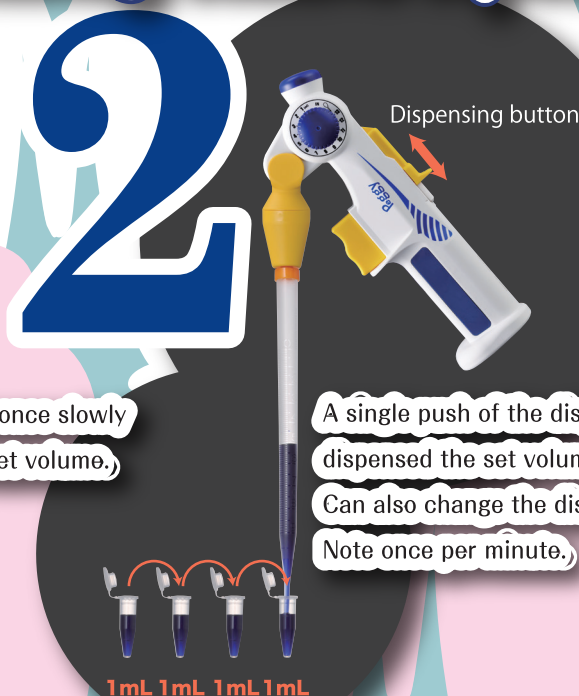

A single push of the dispensing button
dispensed the set volume.
Can also change the dispensing capacity
Note once per minute.

Setting aspiration volume



Aspiration can be set in every
1mL for setting range 1-13mL.


Note: 13mL can be dispensed in case that
Watson 10mL Serological Pipette is used.

Setting dispensing volume




Dispensing can be set in every
0.5mL for setting range
0.5-5mL.

Flexible nozzle-angle

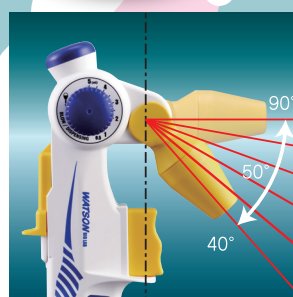

Adjustable nozzle angle in
every 10 degrees up 50 degrees.

Simple maintenance

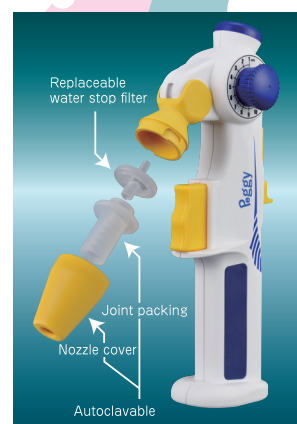

Easy to remove & set spare parts
- Nozzle cover: Autoclavable
- Joint packing: Autoclavable
- water stop filter: Replaceable

Precaution in case of using 5mL serological pipette:

Please DO NOT set aspiration volume exceeding 5mL
in case that 5mL serological pipette is used.
It may cause an accidental liquid suction leading to
malfunction on Peggy.

Cat.No.	Item	Aspiration Volume		Discharge Volume		Packaging
		max.	min.	max.	min.	
PG-01	PEGGY-01 Repetitive Pipette "Peggy"	13 mL	1 mL	5 mL	0.5 mL	1 unit / box

Something Different.
WATSON® BIO LAB
MADE IN JAPAN SINCE 1988

<https://watsonbiolab.com>

Sales agency:
WATSON CO., LTD.

Export Division:
2-2-7, Murotani Nishi-ku Kobe, 651-2241, JAPAN
TEL 81-78-991-4489 FAX 81-78-991-4491

Head Office:
6-57-12, Minamisenju Arakawa-ku, Tokyo,
116-0003, JAPAN
TEL 81-3-5615-3591 FAX 81-3-5615-3592

E-mail:
tcr@watson.co.jp

Manufacturer:
FUKAEKASEI co., LTD.

※The specification and the price etc. of the product might change without a previous notice.